

# How to set up data sharing (for staff)

- Log into a beamline computer (oswin, grimlok, etc) and run the following commands
- `source /home/dm/etc/dm.setup.csh`
- `dm-27id-daq --data-directory=/home/beams/RIXS/Data/2019-1/Slot1-commissioning --experiment=2019Feb12 --duration=168h --esaf-id=202207`

Directory where spec is writing

Experiment name on Globus. Strongly recommend naming by starting date to ensure unique name.

Continues updating for 1 week (168 hours)

Experiment ID from ESAF form. This gives reading permission to users listed on the ESAF

## APS ESAF - Experiment Hazard Control Plan Report

Printed date: 02/13/2019

**PEN:** 27-IDB-2019-164488  
**ID Start Date:** 02/12/2019 08:00 AM  
**Spokesperson:** Bertinshaw  
**Title:** Magnetic excitations in hole doped Sr2Ir[1-x]Rh[x]O4: Links to cuprate superconductors

**Experiment ID:** 202207 (GUP)  
**ID End Date:** 02/18/2019 08:00 AM  
**GUP ID:** 59179